

ACINAR CELLS



ISOLATION OF ADIPOCYTES OR ADIPOSE DERIVED STEM CELLS WITH COLLAGENASE NB

In the following, application details for isolation of adipocytes from human or murine adipose tissue are presented.

Species	Enzyme	Conditions	
Human	Collagenase NB 4	0.2-0.3 PZ U/ml*, 30-60 min, 37 °C	
Murine	Collagenase NB 4	0.2-0.3 PZ U/ml**, 30-60 min, 37 °C	

Clinical Applications**

Species	Enzyme	Conditions	
Human	Collagenase NB 6	0.2-0.3 PZ U/ml*, 30-60 min, 37 °C	

*0.1-0.15 PZ U/ml final concentration after addition of adipose tissue

**For further information on GMP grade Collagenase NB please [contact us!](#)

[Collagenase NB 4 Standard Grade \(cat. no. S1745401\)](#) has a collagenase activity of ≥ 0.1 PZ U/mg and a balanced mix of proteolytic side activities.

[Collagenase NB 6 GMP Grade \(cat. no. N0002779\)](#) has a collagenase activity of ≥ 0.1 PZ U/mg and a balanced mix of proteolytic side activities. It is manufactured according to GMP guidelines and sterile according to European Pharmacopoeia.

Isolation conditions

The suggested Nordmark Collagenase NB concentrations and isolation conditions should be regarded as starting points. Collagenase NB dosage is recommended in PZ units to ensure consistent results.

For further information on optimization of tissue digestion please contact us at (631) 348-0333 or collagenase@crescentchemical.com or visit our website at www.crescentchemical.com.

Ordering:

Crescent Chemical Co., Inc.
2 Oval Dr.
Islandia, N.Y. 11749 U.S.A.
Phone (631) 348-0333 Fax: (631) 348-0913
E-Mail: collagenase@crescentchemical.com
www.crescentchemical.com